

# MC-Media Pad™ “*Salmonella*” instruction manual

## Easy and accurate dry culture system for Microbial Counts

### ◆ BACKGROUND

For hygiene control, it is important to determine the microbial count in foodstuffs and the process environment. MC-Media Pad *Salmonella* is intended to detect *Salmonella* using a special medium composition and specific chromogenic substrate for this pathogen. MC-Media Pad pre-sterilized, ready-to-use dry culture devices simplify testing and minimize the quantity of waste. MC-Media Pad is composed of a unique adhesive sheet, a test pad coated with medium and water absorption polymer, and a transparent cover film. MC-Media Pad is made by ISO 9001 certified factory.

### ◆ TEST PRINCIPLES

MC-Media Pad *Salmonella* is coated with a selective medium and a chromogenic substrate for specific determination. Once the liquid sample is inoculated onto the test pad, the sample diffuses through the whole pad by capillary action. The medium re-constitutes automatically. If *Salmonella* is present, it grows as light-blue colored colonies on the test pad.

### ◆ CONTENTS and STORAGE

- 1000 plates ... code **SK06A25** (25 plates x 40)  
**SK06A10** (10 plates x 100)
- 100 plates ... code **SK06B25** (25 plates x 4)  
**SK06B10** (10 plates x 10)

This kit should be stored between 2-15°C. (Refrigerated)

### ◆ MATERIALS REQUIRED BUT NOT PROVIDED

- Incubator (35 ± 1 °C)
- Stomacher or Blender
- Sampling bag (Recommended for Stomacher; bag with filter to eliminate food debris)
- Pipette or Pipettor and pipette tips
- Butterfield's Phosphate Buffer, Phosphate Buffered Saline, Saline or appropriate diluents
- Pre-enrichment (e.g. BPW) or selective enrichment (e.g. RV broth, TT broth)

### ◆ SAMPLE PREPARATION

- **For solid food stuffs**  
Homogenize a 10-g test portion in 90 mL of Butterfield's Phosphate Buffer, Phosphate Buffered Saline, Saline or appropriate diluents with a stomacher. If necessary, make a 10-fold serial dilution.
- **For water, liquid food stuffs, or swab test sample**  
Sample can be applied directly or diluted with Butterfield's Phosphate Buffer or appropriate diluents as for solid foodstuffs. If necessary, pH of sample should be adjusted to neutral (pH 7.0±0.2).

### ◆ TEST PROCEDURE

- **General Operation**
  1. Open the aluminum bag, and remove the MC-Media Pad.  
If necessary, write information on the cover film.
  2. Lift the cover film and drop 1.0 mL of sample solution onto test pad.
  3. Replace the cover film and lightly press the edges of film to seal. (It is recommended to lift the cover film diagonally for easy and sure re-sealing.)
  4. Incubate test plate at 35 ± 1°C for 24 ± 2 hours.
- **Other Application**  
MC-Media Pad is also available for Wiping/Stamping technique, Membrane filter method, and Airborne falling bacteria test. MC-Media Pad website provides detailed information.  
(<https://www.jnc-corp.co.jp/MC-MP/>)

### ◆ INTERPRETATION

Count all light-blue colored colonies as *Salmonella* regardless of strength of color. Purple or red colored colonies is not *Salmonella*. For large number of colonies, colony counts can be estimated by counting colonies in one grid square and multiplying by 20.

If more than 10<sup>4</sup> microbes are grown, the entire test pad may appear as stained, and it may appear that no individual colonies were formed. If this is the case, dilute the sample further and re-test. If necessary, a target colony can be picked up with a sterile needle from the test pad for further analysis.

### ◆ PRECAUTIONS

1. The test is designed for use by quality control personnel and others familiar with testing samples potentially contaminated with *Salmonella*.
2. Read this instruction manual carefully before use.
3. After opening the aluminum bag, unused plates should be stored in the aluminum bag sealed with tape, and kept in a cool (2-15°C) environment. After opening, use all plates within 1 month.
4. Do not expose unused plates to sunlight or ultraviolet light.
5. Do not use a discolored or damaged plate.
6. A wrinkle on the test pad should not affect detection.
7. Small fragments of fabric on or around the test pad should not affect detection.
8. Do not use the plates after the expiration date. The quality of an expired plate is not warranted.
9. The measurement range is less than 300 cfu/plate. If more than 300 cfu/plate are counted, further dilution is recommended.
10. The presumptive *Salmonella* detected on MC-Media Pad “*Salmonella*” should be confirmed to be *Salmonella* by serological test or other tests.
11. In case of applying enzyme-containing foods (e.g. raw foods or liver), the entire test pad may appear as stained.
12. The nature (high viscosity food or food dye) of food may affect test usage or results. In that case, the causes need to be eliminated by dilution or other means.
13. The used kit must be sterilized by autoclaving or boiling, and then disposed according to local regulations for waste.

### ◆ LIMITATION of WARRANTY

The Products are covered by the applicable JNC Corporation standard warranty. NO OTHER EXPRESS OR IMPLIED WARRANTY IS MADE WITH RESPECT TO THE PRODUCTS. JNC EXPRESSLY EXCLUDES THE IMPLIED WARRANTIES OF MERCHANTABILITY AND OF FITNESS FOR A PARTICULAR PURPOSE. If product is defective, JNC and JNC's authorized distributor will provide a replacement or refund at the purchase price.

### ◆ CONTACT and FURTHER INFORMATION

#### JNC Corporation Life Chemical Division

2-2-1 Otemachi, Chiyoda-ku, Tokyo 100-8105 Japan  
TEL:+81-3-3243-6225,FAX:+81-3-3243-6219  
E-mail: [mcmp@jnc-corp.co.jp](mailto:mcmp@jnc-corp.co.jp)

Manufactured by **JNC CORPORATION**